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Electronic components must not be disposed of in the domestic waste at the end of their service life. Used electronic devices contain harmful substances that can cause damage to the environment or human health. End users are legally obliged to take used electric and electronic devices to a licensed collection point.

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Trademark index:

VACUU·LAN® (US-Reg.No 3,704,401), VACUU·BUS®, VACUU·CONTROL®, VACUU·SELECT® (US-Reg.No 5,522,260), VARIO® (US-Reg.No 3,833,788), VACUUBRAND® (US-Reg.No 3,733,388), VACUU·VIEW®, GREEN VAC® (US-Reg.No. 4,924,553), VACUU·PURE® (US-Reg No. 5,559,614) and also the shown company logos are registered trademarks of VACUUBRAND GMBH + CO KG in Germany and/or other countries.

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Safety information!

NOTICE

The coolant valves 20676010 and 20676013 are to be used only in combination with the vacuum controllers CVC 2II, CVC 2000, or VNC 1, or with the Vacuum Management System VMS Module A. The coolant valve 20674220 is designed to be connected to the VACUUBRAND BUS system VACUU•BUS.

Remove all packing material from the packing box. Remove the product from its packing-box and retain all packaging for inspection until the equipment is inspected and tested. Inspect the equipment promptly and carefully. If the equipment is damaged, notify the supplier and the carrier in writing within three days. State the item number of the product together with the order number and the supplier's invoice number.

Use the equipment **for the intended use only**, i.e. to control the coolant flow at the inlet of the cooling system.

Comply with all applicable and relevant safety requirements (regulations and guidelines). Implement the required actions and adopt suitable safety measures.

Ensure that **maintenance** is done only by suitably trained and supervised technicians. As laid down in the statutory regulations (occupational, health and safety regulations and regulations for environmental protection), equipment returned to the manufacturer can be repaired only under certain conditions.


WARNING

- ☞ **Do not use the equipment if it is damaged.**
- ☞ Use only genuine spare parts and accessories. Otherwise safety and performance of the equipment as well as the electromagnetic compatibility of the equipment might be reduced.
- ☞ Pay attention to the permissible maximum ambient temperatures.
- ☞ Check that the system is mechanically stable and that all fittings are secure.

CAUTION

- Check that the power source and the equipment match in voltage, phase, and frequency (see rating plate).
- Ensure that the coolant is flowing in the correct direction (see arrow on valve housing).
- Ensure that the materials of the wetted parts are compatible with the media, see "Technical data".

Coolant valve 230 V~ 50/60 Hz, IEC plug	20676010
Coolant valve 230 V~ 50/60 Hz, CEE plug	20676014
Coolant valve 230 V~ 50/60 Hz, UK plug	20676012
Coolant valve 24 V= VKW	20676013
Coolant valve 24 V= VKW-B	20674220

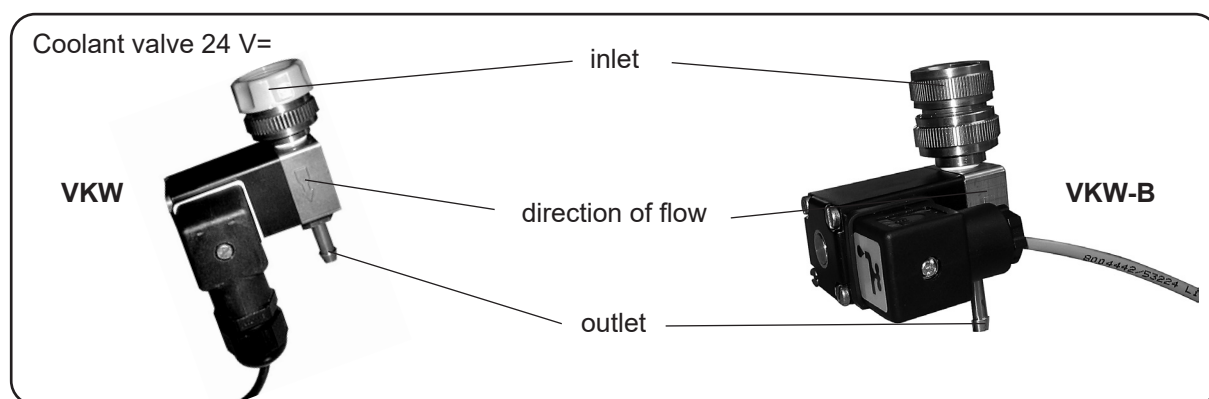Attention! Additional notes concerning customized equipment:

Material No. 22610867: hoses nozzles (DN 6) at water inlet and water outlet.

Material No. 22611381 / 22613924: threaded connections 1/2" (3/4") at water inlet and water outlet

Technical data

Type	20676010 20676012 20676014	20676013 20674220
Maximum permissible range of supply voltage	230 V~ ($\pm 10\%$) 50/60 Hz	24 VDC $\pm 10\%$
Power draw (operation)	8 W	2 W
Electrical connection	20676010: IEC plug (VMS Module A) 20676012: UK plug 20676014: CEE plug	20676013: male connector 5-pole* 20674220: connector for VACUUBRAND BUS system VACUU•BUS
Rated duty	continuous operation	
Valve currentless	valve closed	
Nominal diameter	2.0 mm	1.5 mm
Throughput (Kv value water**)	2.0 l/min	1.3 l/min
Max. permissible water pressure	10 bar	5 bar
Inlet connection	G3/4" with reducing piece G1/2"	
Outlet connection	G1/8" with hose nozzle DN 6 mm	
Permissible ambient temperature	0°C to +40°C	
Permissible media temperature	-10°C to +40°C	
Dimensions L x W x H (without cable)	92 mm x 73 mm x 72 mm	100 mm x 65 mm x 72 mm
Cable length	2.5 m	2 m
Weight	0.57 kg	0.45 kg

* cable with integrated Y-socket-plug connector to connect a coolant valve and an in-line valve to the controllers CVC 2[®], CVC 2000, or VNC 1 or to the Vacuum Management System VMS Module A.

** measured at +20°C, with 1 bar water pressure at the inlet and free water flow at the outlet

Components	Wetted materials
Water connection	brass
Hose nozzle	brass
Reducing piece	brass
Seal ring R1/8"	PVC
Flat packing R1/2"	NBR
Flat packing R3/4"	FPM

We reserve the right for technical modifications without prior notice!

Spare parts

1x Flat packing R3/4"	20610603
1x Flat packing R1/2"	20610604

Documents are only to be used and distributed completely and unchanged. It is strictly the users' responsibility to check carefully the validity of this document with respect to his product.

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DECLARATION OF CONFORMITY – China RoHS 2

VACUUBRAND GMBH + CO KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a “Product Conformity Assessment” (PCA) procedure was performed. As defined in GB/T 26572 the “Maximum Concentration Value” limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(+VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmentally Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user’s assets. The Environmentally Friendly Use Period for VACUUBRAND products is 40 years.

MATERIAL CONTENT DECLARATION FOR VACUUBRAND PRODUCTS						
有毒有害物质或元素 Hazardous substances						
部件名称 Part name	铅 Pb	汞 Hg	镉 Cd	六价铬 Cr(+VI)	多溴联苯 PBB	多溴二苯醚 PBDE
包装 Packaging	0	0	0	0	0	0
塑料外壳 / 组件 Plastic housing / parts	0	0	0	0	0	0
真空油 Vacuum oil	0	0	0	0	0	0
电池 Battery	0	0	0	0	0	0
玻璃 Glass	X	0	0	0	0	0
电子电气组件 Electrical and electronic parts	X	0	0	0	0	0
控制器 / 测量设备 Controller / measuring device	X	0	0	0	0	0
金属外壳 / 组件 Metal housing / parts	X	0	0	0	0	0
电机 Motor	X	0	0	0	0	0
配件 Accessories	X	0	0	0	0	0
此表格是按照SJ/T 11364-2014中规定所制定的。 This table is created according to SJ/T 11364-2014.						

VACUUBRAND®

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
 O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
 X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Batteries, glassware and accessories might not be content of the enclosed device and may have its own EFUP-marking and/or might be maintaining parts with changing EFUP-marking.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (Cr+VI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by VACUUBRAND may enter into further devices (e.g., rotary evaporator) or can be used together with other appliances (e.g., usage as booster pumps).

With these products and appliances in particular, please note the EFUP labeled on these products.

VACUUBRAND will not take responsibility for the EFUP of those products and appliances.

Place, date: Wertheim, 06 September 2022

(Dr. Constantin Schöler)
 Managing Director

ppa.

(Jens Kaibel)

Technical Director

VACUUBRAND GMBH + CO KG



Alfred-Zippe-Str. 4
 97877 Wertheim
 Germany

Tel.: +49 9342 808-0

Fax: +49 9342 808-5555

E-Mail: info@vacuubrand.com

Web: www.vacuubrand.com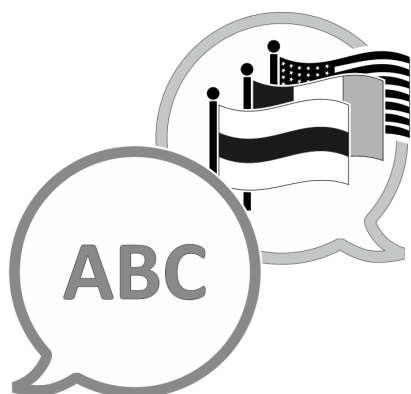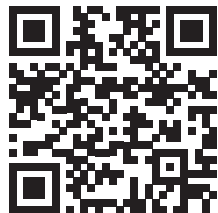

[VACUUBRAND > Support > Manuals](#)

Manufacturer:

VACUUBRAND GMBH + CO KG
Alfred-Zippe-Str. 4
97877 Wertheim
GERMANY

Phone:

Head office +49 9342 808-0
Sales +49 9342 808-5550
Service +49 9342 808-5660

Fax: +49 9342 808-5555

Email: info@vacuubrand.com

Web: www.vacuubrand.com