AUSBAUSÄTZE UPGRADE KITS

20699979 AK

20699948 EK

Montageanleitung / Mounting instructions

Unbedingt beachten!

VORSICHT

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Safety information!

CAUTION

- Ensure that installation work is done only by a suitable trained and supervised technician. Obey local and national safety requirements. Ensure that the technician is familiar with the safety procedures which relate to the substances processed by the pumping system. Ensure that the device is decontaminated before repair and that you take adequate precautions to protect people from the effects of dangerous substances if contamination has occurred.

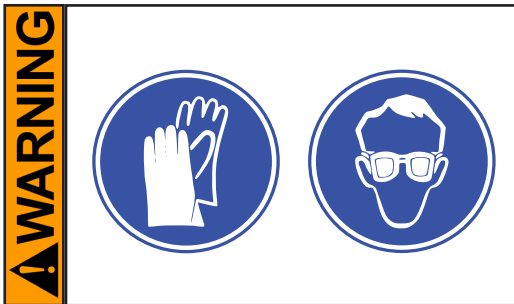
DANGER


- ➔ **Never operate the pump if covers or other parts of the pump are disassembled.**
- ➔ Before starting your work, **disconnect the electrical power cord. Wait two minutes** after isolating the equipment from AC power to allow the capacitors to discharge.
- ➔ Ensure that the pump cannot be operated accidentally.
- ➔ **Note:** The pump might be contaminated with the process chemicals that have been pumped during operation. Ensure that the pump is decontaminated before starting any installation work.
- ➔ Avoid the release of pollutants.

WARNING

- ➔ **Never operate a defective or damaged pump.**
Do not reuse damaged parts.
- ➔ Use only **original manufacturer's spare parts and accessories.** Otherwise the safety and performance of the equipment, as well as the electromagnetic compatibility of the equipment might be reduced.
- ➔ Read the **instructions for use** for the equipment carefully, especially sections "Safety informations" and "Use and operation". If necessary order "Instructions for use" from VACUUBRAND or download at www.vacuubrand.com.

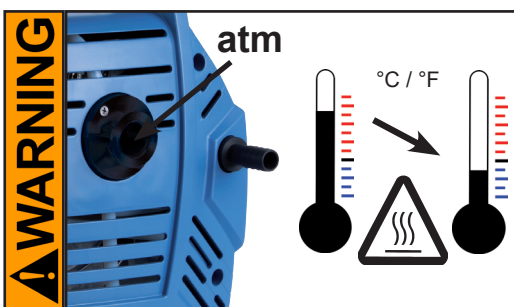- ➔ Take adequate precautions to protect people from the effects of dangerous substances that may have contaminated the pump and may be released upon disassembly. Use appropriate protective clothing, safety goggles and protective gloves.
- ➔ Comply with applicable regulations when disposing of chemicals. Take into consideration that chemicals may be contaminated.

- ➔ Allow sufficient cooling of the pump **before starting installation work.**
Vent the pump and isolate it from the vacuum system before you start your work. Drain condensate, if applicable.

Before starting installation work check that the required tools and parts are available and of the correct type.

Check the operating sequence mentally on feasibility, safety requirements and consequences on safety and function of the equipment.

The pictures may show other versions of pumps. This does not change the method of assembling the separator or the emission condenser kit.

CAUTION

- **Check operability and safety after installation work has been completed.**
In case of problems return the device to the factory for inspection or repair if necessary.
- Pay attention to the safety symbol "hot surfaces" on the equipment. Depending on operation conditions and ambient conditions dangers due to hot surfaces may arise. Hot parts may cause burns if touched. Adopt suitable measures to prevent any danger arising from hot surfaces. Ensure that hot surfaces of the pump do not cause burns. Provide a suitable contact guard if necessary.
The surface temperature of the glass components at the exhaust waste vapor condenser in particular might be elevated in case of permanent high gas flow. The temperatures that are produced during operation may cause burns. Avoid direct contact with the surface. Provide a suitable contact guard if necessary.

In order to comply with laws (occupational, health and safety regulations, safety at work law and regulations for environmental protection) vacuum pumps, components and measuring instruments can only be returned when certain procedures (see section "**Notes on return to the factory**" in the "Instructions for use") are followed.

Scrapping and waste disposal

Dispose of the pump and any components removed from it safely in accordance with all local and national safety and environmental requirements.

Particular care must be taken with components and waste oil which have been contaminated with dangerous substances from the process.

 You may authorize us to dispose the equipment **at your expense**.

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